# MC 1 MC 2 – MC 3 MC 4

## Manual Capsule Filling Machines

MACHINE	BATCH Q.TY	PRODUCTION / H	CAPSULE SIZES	ACCESSORIES	DIMENSION s LxWxH CM	WEIGHT KG.
MC 1	50	1200	000/5	PRESSORS,	19x16x7	1
MC 2	100	1200	000/5	HOPPERS,	18x22x10	2
MC 3	60	1500	00/4	VIBRATING PLATE	19x18x7	1
MC 4	120	1500	00/4		18x25x10	2

MACHINE	POWER CONSUMPTION KW	POWER SUPPLY KW	DIMENSIONS LxWxH CM	WEIGHT KG.
PV1	0,07	110/230 SINGLE PHASE	24X17X10	3

### Macchine manuali per capsule

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### Manual Capsule Filling Machines

The manual capsule orientation device MC 1, capable of processing capsules from size 000 to 5, can align 100 capsules in two steps. In turn, the manual filling machine MC 2 (also capable of processing capsule sizes from 000 to 5) can open, fill and close up to 100 capsules at a time. The capsule opening system has been tried and tested for more than twenty years and its building materials have been constantly updated, including recently: The kit to switch capsule size is made of a special patented resin (ERTACETAL), while stainless steel is used for the parts in contact with the product. The frame is made of anodized hard aluminum, to facilitate cleaning.

It is possible to switch size using the transformation kit.

The set of orientation/filling machines MC 3 MC 4 differs only in the total number of capsules, 120 instead of 100.

Accessories are available for purchase to facilitate the preparation: both fixed and adjustable powder hoppers, pressure pushers in different sizes, and vibrating table PV 1 to facilitate the smooth and even descent of the powders into the capsules.

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